



AWRI
COMMERCIAL SERVICES
SO MUCH MORE THAN A GREAT LAB

Fact Sheet

Sequencing yeast in natural ferments



Sequencing service

AWRI Commercial Services can apply next-generation sequencing technologies to provide information about the dominant yeast species present in 'natural' ferments.

Why is it important?

Natural, spontaneous, wild and uninoculated are several terms used to describe a wine fermentation that has not been inoculated with a commercial yeast strain. The population of yeast species is dynamic throughout an uninoculated fermentation due to the range of initial microflora present and changing conditions as the fermentation progresses. The yeast population is most diverse for the first few days of fermentation, after which *S. cerevisiae* tends to dominate as alcohol levels increase. This service will provide insight into the microflora present

during fermentation that may contribute characteristic flavours to wine.

How does it work?

The genomic DNA is extracted from a ferment sample provided by the winery. Certain sections of the DNA are amplified and then sequenced which then provides information about the species present.

What does it involve?

The AWRI will provide a sampling kit containing a small esky, ice pack and sample tubes. The client should take a 50 mL sample of the ferment and then allow it to cold settle in the fridge overnight. The next day the liquid should be poured off and the solids that have settled out left in the tube and frozen. The client should then send the esky containing the sample and ice pack to the AWRI by



AWRI



AWRI
COMMERCIAL SERVICES
SO MUCH MORE THAN A GREAT LAB

Fact Sheet

courier. It is recommended that samples are taken during early stages of fermentation, for example at crushing and then after one, two and three days of fermentation. However a single sample in the early stages of fermentation is still able to provide a snapshot of species present.

What will the results show?

Results from the analysis will provide a list of the yeast species present in the sample and approximate proportions of different species.

How much does it cost?

Price depends on the number of samples being sequenced. Please contact us for an individualised quote.

Additional services available

Genetic characterisation of natural ferments to the strain level

In this service, all the genetic material is sequenced to the strain level. Such analysis can determine if a ferment contains commercial yeast strains and will also characterise strains that may be unique to a particular winery.

Isolation of dominant yeasts and cryogenic storage of them in the [AWRI Wine Microorganism Culture Collection](#)

This may act as insurance for future vintages, preserving microorganisms that are likely to have an impact on wine characters.

Contact

For further information, please contact:

Commercial Services

Phone 08 8313 6600 Fax 08 8313 6601

Email CommercialServices@awri.com.au

Address Wine Innovation Central Building,
Corner of Hartley Grove & Paratoo Rd,
Urrbrae (Adelaide), SA 5064